SCIENTIFIC EQUIPMENT

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Description: Ash Fusibility Determinator is used to determine the behavior of Coal and Coke ash at temperature 900 – 1500 °C. Design of most coal combustion and coal conversion equipment anticipates that the ash either remains. The result can be report of IT or DI (Initial Deformation Temperature), ST (Spherical Temperature), HT (Hemispherical Temperature), and FT (Fluid Temperature).

Sample types: Coal and Coke ash

SURFACE AREA AND MICROPORE ANALYZER (BET)

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Description: Analyzer for surface area and pore size distribution of particles automatically, based on static the measurement of gas volume which is adsorbed on the surface of particles with static volumetric gas adsorption method and can measure pore size distribution in the range of 3.5 angstroms – 5000 angstroms. It has three pressure transducer systems which are Transducers 1000 mmHg with 0.001 mmHg resolution, Transducers 10 mmHg with 0.00001 mmHg resolution, Transducer 0.1 mmHg with 0.0000001 mmHg resolutions. Reports include summary report, isotherm report, single and multipoint BET Surface area, Langmuir Surface area,

total pore volume, BJH adsorption and desorption, Dollimore-Heal adsorption and desorption, t-Plot method for micropore volume and area, Sample log, Horvath-Kawazoe, Dubinin, DFT, NLDFT, MP reports.

Sample types: It is ideally suited for the characterization of pharmaceuticals, ceramics, activated carbons, carbon black, catalyst, paints and coatings, projectile propellant, medical implants, electronics, cosmetics, aerospace, geoscience, nanotubes, and fuel cells, *etc.*

HIGH THROUGHPUT SURFACE AREA AND POROSITY ANALYZER (BET)

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Description: Surface Area and Porosimetry Analyzer is Static Volumetric Gas Adsorption Method. This instrument can measure specific surface area minimum $0.01 \text{ m}^2/\text{g}$ by Nitrogen gas, pore volume not less than $4 \times 10^{-6} \text{ cm}^3/\text{g}$. There are 6 adsorptive gas inputs.

Sample types: It is ideally suited for the characterization of Pharmaceuticals, Ceramics, Activated Carbons, Carbon Black, Catalyst, Paints and Coatings, Projectile Propellant, Medical Implants, Electronics, Cosmetics, Aerospace, Geoscience, Nanotubes, and fuel cells, etc.

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Description: DMA is used for analyzing the physical properties of materials that are related to temperature and time such as Glass Transition, Temperature, Storage Modulus, Loss Modulus and Tan Delta. The analysis can be made through the methods of Dynamic, Tension, Bending, Compression, and Shear in the range of -190 °C to 600 °C.

Sample types: DMA can analyze the Samples which are plastic, polymers, rubber and film, etc.

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Description: Rheometric Scientific DMA is used for analyzing the physical property of materials that are related to temperature and time such as Glass Transition, Temperature, Storage Modulus, Loss Modulus and Tan Delta. The analysis can be made through the methods of Dynamic, Tension, Bending, Compression, and Shear in the range of -150 to 500 °C The innovative, material pockets allow powdered or non-self-supporting materials, such as powdered drugs, gels, natural products like tea, coffee, herbs, etc.

Sample types: plastic, polymers, rubber, film, powder, gels, and natural products such as tea, coffee, herbs, etc.

DIFFERENTIAL SCANNING CALORIMETER (DSC1)

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Description: DSC is used for analysis the physical property of materials such as boiling point, melting point which it's changing temperature point ranges from 20°C - 500°C . Only using 2 mg - 10 mg of sample for testing.

Sample types: DSC7 can analyze plastic, polymer, drugs, animal protein, and clay, etc.

DIFFERENTIAL SCANNING CALORIMETER (DSC2)

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Description: DSC is used for analysis the physical property of materials such as boiling point, melting point which its changing temperature point ranges from $-170\text{ }^{\circ}\text{C}$ to $500\text{ }^{\circ}\text{C}$. Only using 2 mg - 10 mg of sample for testing.

Sample types: The equipment can analyze plastic, polymer, drugs, animal protein, and clay, etc.

DIFFERENTIAL SCANNING CALORIMETER (DSC3)

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- Melting behavior
- Crystallization and nucleation
- Polymorphism
- Liquid-crystalline transitions
- Phase diagrams and composition
- Glass transitions
- Reactivity
- Reaction kinetics
- Curing
- Stability
- Miscibility
 - Effects of plasticizers
- Thermal history
- Heat capacity and heat capacity changes
- Reaction and transition enthalpies
- Purity

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Description: DSC is used for measures the enthalpies associated with transitions and reactions and the temperatures at which these processes occur. The method is used for the identification and characterization of materials. Differential scanning calorimetry (DSC) is fast and very sensitive. Sample preparation is easy and requires only small amounts of material. The technique is ideal for quality control, material development and material research. DSC is the method of choice to determine thermal quantities, study thermal processes, and characterize or just simply compare materials. It yields valuable information relating to processing and application conditions, quality defects, identification, stability, reactivity, chemical safety and the purity of materials. The method is used to analyze and study materials such as thermo-plastics, thermosets, elastomers, composite materials, adhesives, foodstuffs, pharmaceuticals and chemicals. Thermal events and processes that DSC can be determined is

- Melting behavior
- Crystallization and nucleation
- Polymorphism
- Liquid-crystalline transitions
- Phase diagrams and composition
- Glass transitions
- Reactivity
- Reaction kinetics
- Curing
- Stability
- Miscibility
- Effects of plasticizers
- Thermal history
- Heat capacity and heat capacity changes
- Reaction and transition enthalpies
- Purity

Sample types: Used to analyze material samples such as Polymer (Thermoset/Thermoplastic), flim, rubber, pharmaceutical and chemical, *etc.* and biological samples such as Protein product (Fish, Egg), *etc.*

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Description: DTA is used for analyzing the physical property of materials such as boiling point, melting point which it's changing temperature point ranges from $50\text{ }^{\circ}\text{C}$ – $1300\text{ }^{\circ}\text{C}$. Only using 2 mg - 10 mg of sample for testing.

Sample types: DTA can analyze plastic, polymer, drugs, animal protein, and clay, *etc.*

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Description: It is used for measuring the particle size and the distribution of the samples which are powder and suspension. It can analyze in both the liquid and air medium. It can measure the particle size from 0.04 micrometer to 2,000 micrometer in liquid medium and 0.4 micrometer to 2,000 micrometer in air medium.

Sample types: The samples are powder, liquid and suspension forms such as concentrated latex, synthetic latex and chemicals in industrial rubber gloves.

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Description: It is used for measuring the particle size and the distribution of the samples which are powder and suspension. It can analyze in both the liquid and air medium. It can measure the particle size from 0.01 micrometer to 2,000 micrometer in liquid medium and 0.1 micrometer to 2,000 micrometer in air medium.

Sample types: The Samples are powder, liquid and suspension forms such as concentrated latex, synthetic latex and chemical in industrial rubber gloves.

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Description: STA8000 is used for analyzing the physical property of materials such as boiling point, melting point and detecting the missing weight volume because of the disintegration of composites in the Samples when temperature changes. The changing temperature point of STA8000 ranges from 25°C to 1300 °C. Only using 2 mg - 10 mg of sample for testing.

Sample types: The materials studied include plastic, polymer, drugs, animal protein, clay, soil, rubber, wood, ceramic and etc.

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Description: It is used for detecting the missing weight volume because of the disintegration of composites in the samples when temperature changes in the range of 50 °C to 1300 °C.

Sample types: The materials studied including of soil, rubber, wood, ceramic, plastic, and etc.

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Description: It is real-time measurement and analysis of sample weight change while subjecting it to a controlled continuously of the optimization in temperature time and atmosphere. An application is characterizing inorganics, analyzing polymers mineral or coal. Suitable for liquid, solid, powder, film, and fiber etc. Temperature range of a single furnace start from ambient to 1,000 °C with embed heater in furnace wall. The ramp rate of heating is up to 50°C/min. The instrument consist of the balance readability to 0.0001 g and a standard weight-loss for an inert Sample with a precision of $\leq 0.02\%$ standard deviation.

Sample types: The samples can be both solids and liquids such as rice husk, oil, palm shell, coal or biomass product, *and etc.*

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Description: TCA is used for analyzing thermal conductivity thermal diffusivity and specific heat of many kinds of samples.

Sample types: The samples can be solid, liquid, film, powder, cream and etc. Samples must be smooth and homogeneous with size at least 5 cm x 5 cm x 1 cm. For film sample and paper size are 50 mm x 50 mm x 0.02 mm and liquid Sample volume at least is 100 mL.

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Description : Zeta is used to detect the zeta potential values of colloid with zeta potential values between -150 mV and $+150\text{ mV}$ as well as measure the size of nanoparticles ranging between 8.4 nanometer and 30 microns.

Sample types: The samples tested can be aqueous (water, methanol, and ethanol) and non-aqueous: Samples can be powder, liquids, or colloid such as pigment, ceramics, soil, pharmaceutical products, and etc.

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Description: It is used for rheological measurements with control the temperature such as viscosity, elastic properties, deformation of fluid and it can analyze flow behavior of many types of fluid.

Sample types: Viscous fluid such as oil, cream and gel, etc.

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Description: It is used to measure the volume, density and specific gravity of the sample, by using the gas displacement method to measure volume and calculate density. Inert gases, such as helium or nitrogen are used as the displacement medium. The instrument provides high-speed, high-precision volume measurements and true density calculations.

Sample types: The samples can be a wide varieties of powders, solids and liquid such as animal feeds, cements, ashes, *and etc.*

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Description: Density analyzer is used to measure the actually density of the sample based on the electromagnetically induced oscillation of a U-form tube made of glass. A magnet is fixed to the U-tube and a transmitter induces the oscillation. The period of oscillation is measured by a sensor.

Sample types: The samples must be dissolved in organic solvents and deionized water, except for the viscous samples which get clump easily, such as latex, honey and gel, etc.

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Description: The equipment measures both reflected and transmitted color as well as transmission haze and meets CIE, ASTM and USP. The results of this instrument are absorbance, brightness, color, haze, K/S, metamerism, opacity, reflectance, strength, transmittance, whiteness, yellowness, and gray scale, etc.

Sample types: This instrument can measure the color of samples in both solid and liquid forms from opaque to translucent materials, such as film, oil powder, waste water, ceramics, rubber, and natural materials, etc.

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Description : The analyser measures the permeation rate in chemical safety natural and synthetic rubber gloves. The chemicals are detected using a built-in FID. The analyser measures the breakthrough of the test chemicals in permeation measurement cells according to DIN EN 374-3.

Sample types: Natural and synthetic rubber gloves.

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Description: pH meter is a device that can measure acidity - alkalinity in the solution. pH can be measured from 0.00 to 14.00. The pH reading is 0.01 pH.

Sample types: Liquid Samples such as drinking water, raw water, and wastewater, etc. Solid Samples must be prepared in to solution before testing.

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Description: Electrical conductivity can be measured by using the resistance of the solution between the anode and the cathode. Which when dipping both types of probes into solution, negative ions will move towards the anode. And the cation will move towards the cathode. The movement of ions results in a rising of current. The current occurs more or less, depends on the resistance of the solution between two electrodes.

Sample types: Liquid and solid samples such as drinking water, raw water, wastewater, soil, meat and herbs, etc.. Solid Samples must be prepared in to solution before testing.

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Description : Viscometer is an instrument that used to measure a viscosity of a liquid with using the spindle rotating in the fluid. It can read the viscosity of a liquid in centipoise (cP) unit.

Sample type : Liquid samples such as latex, honey, and milk, etc.

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Description: It used to measure a viscosity at 40°C according to ASTM D 445. Kinematic Viscosity of liquid samples was measured by measuring the time of the liquid flowing under the gravitation through the viscometer tube. The viscosity in the sample is the result of the time that the fluid flows from the upper mark to the lower mark and the viscometer tube constant that used in the sample test.

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Description: Hydrometer is used for testing the specific gravity and the API gravity at 60 ° F and density at 15 ° C in biodiesel by relying on the liquid float principle.

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(ANALYTICAL THE MOLECULAR MASS
AND ORGANIC COMPOUNDS)

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Description: Capability of separating mixture by turning substance into gas at temperature not more than 300 °C with FID detector of GC. The samples analyzed are such as alcohol, pesticide residue, insecticides in crude palm oil or in natural samples. Furthermore, it is capable of qualitative analysis for natural products and fatty acid & methyl ester essential oil. The samples test at temperature not more than 400 °C and they can be volatile and semi volatile substances. The results of the test can be qualitative such as the analysis of the compositions of wood vinegar and essential oil. Quantitative tests are methyl ester in biodiesel oil, essential fatty acid, and essential fatty acid in materials.

Sample type: It can be analyzed including of alcohol and pesticide residues in various samples, natural products, drugs, essential oils, wood vinegar, etc.

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Description: Gas chromatograph mass spectrometer with FID detector is an instrument used to analyze groups of volatile organic compounds and a group of semi-volatile organic compounds using the principle of separation by gas chromatography .The extracted substance can accurately predict the type of components contained in the substance by comparing the mass pattern of the substance of interest with the information contained in the library. It can be used for both quantitative and qualitative analysis.

Sample type : It can be analyzed including of alcohol and pesticide residues in various samples, natural products, drugs, essential oils, and wood vinegar, etc.

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Description: Gas chromatograph Mass spectrometer with CI detector is an instrument used to analyze groups of volatile organic compounds and a group of semi-volatile organic compounds using the principle of separation by gas chromatograph. The extracted substance can accurately predict the type of components contained in the substance by comparing the mass pattern of the substance of interest with the information contained in the library. This can be used for both quantitative and qualitative analysis.

Sample type: Examples that can be analyzed include alcohol and pesticide residues in various samples, natural products, drugs, essential oils and wood vinegar, etc.

GAS CHROMATOGRAPH TRIPLE QUADRUPOLE SPECTROMETER (GC-MS/MS)

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Description : Gas chromatograph Mass spectrometer is a instrument used to analyze groups of volatile organic compounds and a group of semi-volatile organic compounds using the principle of separation by chromatography .The extracted substance can accurately predict the type of components contained in the substance by comparing the mass pattern of the substance of interest with the information contained in the library. It can be used for both quantitative and qualitative analysis.

sample type : alcohol and pesticide residues in various samples, natural products, drugs, essential oils, and wood vinegar, *etc.*

COMPREHENSIVE GAS CHROMATOGRAPH COMBINED WITH TIME-OF-FLIGHT MASS SPECTROMETER (GCxGC-TOF MS)

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Description: The state-of-the-art comprehensive gas chromatograph combined with Time-of-Flight mass spectrometer (TOF-MS) is the powerful tool for separation, analysis, Identification, characterization and fingerprinting of volatile organic compounds in complex samples. The GCxGC system is performed by using two GC separating columns (usually non-polar and polar) connected in series through a modulator. The use of these two separation mechanisms

increases the chromatographic plane. Moreover, the GCxGC system is integrated with automated headspace (HS) and automated headspace solid-phase microextraction (HS-SPME) for a variety of applications. Time-of-Flight mass spectrometer (TOF MS) detector generate full-spectrum data in a very fast up to 500 spectra/s over a mass range of 5 amu - 1000 amu. The spectrum is then compared to known analytes in spectrum libraries. The GCxGC-TOF MS can perform both qualitative and quantitative analyses

Sample type: A sample can be a solid or a liquid such as environmental, petrochemical, foods and beverages, fatty acids, pesticides, plant extracts, natural products, flavour/fragrance, metabolomics and forensic science/toxicology etc.

LIQUID CHROMATOGRAPH-MASS SPECTROMETER / MASS SPECTROMETER (LC-MS/MS)

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Description: LC-MS/ MS instrument is a combination of ultra-high performance liquid chromatograph (UHPLC) with mass spectrometer. UHPLC is used to separate organic compounds from each other and from the complex mixtures or matrices prior to mass spectrometer. Quadruple-time-of-flight mass spectrometer (Q-TOF MS) is a high-resolution mass spectrometer with the ability to measure accurate mass to charge ratios (m/z) with four decimal places. Electrospray ionization (ESI) and atmospheric pressure chemical ionization (APCI) sources are available. The instrument can perform both quantitative and qualitative analyses.

Sample type: solid or a liquid, but the sample must be a solution prior to analysis. LC-MS/MS is widely used in the identification and the characterization of compounds from natural product extracts. Pesticide residues in fruits and vegetables or environmental can be analyzed using LC-MS/MS.

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Description: Gas chromatograph used for analyzing of volatile, VOCs and semi-volatile organic compounds, SVOCs based on the principle of separation between two phases, consisting of stationary phase and mobile phase. The stationary phase is the substances that is coated or packed within the column and mobile phase is carrier gas often use helium gas (He). When the mixtures were injected into the inlet area, which is the area that provides heat is high enough to make the liquid phase change to gas phase and the gas phase of the mixtures will move on into the column by using the carrier gas and the heating of the column with the oven, causing the separation process. Then, the mixture can be separated within the column will be entered into the detector and the result will be displayed in the form of chromatograms.

Sample type: biogas samples consist of hydrogen, nitrogen, methane gas and carbon dioxide gas were kept in a sealed container such as gas sampling bag with a septum can be penetrated through.

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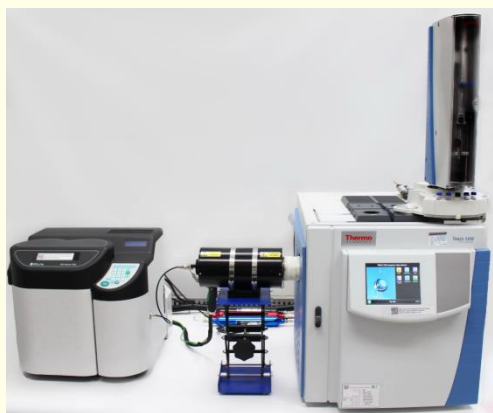
Description: Gas chromatograph is a tool used for analyzing of volatile organic compounds, VOCs and semi-volatile organic compounds, SVOCs based on the principle of separation between two phases, consisting of stationary phase and mobile phase. For the stationary phase of GC technique is the substances that is coated or packed within the column and mobile phase is carrier gas often use is helium gas (He). When the mixtures were injected into the inlet area, which is the area that provides heat is high enough to make the liquid phase change to gas phase and the gas phase of the mixtures will move on into the column by using the carrier gas and the heating of the column with the oven, causing the separation process. Then, the mixture can be separated within the column will be entered into the detector and the result will be displayed in the form of chromatograms.

Sample type : Which can be tested both qualitative and quantitative analysis in many types of sample such as alcohols, organic acids, fatty acids, general oils, biodiesel for FID and organic halide compounds for μ ECD *etc.*

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Description: It uses for analyzing of volatile, VOCs and semi-volatile organic compounds, SVOCs based on the principle of separation between two phases, consisting of stationary phase and mobile phase. When the mixtures were injected into the inlet area, which is the area that provides heat is high enough to make the liquid phase change to gas phase and the gas phase of the mixtures will move on into the column by using the carrier gas and the heating of the column with the oven, causing the separation process. Then, the mixture can be separated within the column will be entered into the detector and the result will be displayed in the form of chromatograms.

Sample type : Which can be tested both qualitative and quantitative analysis in many types of sample such as alcohols, various fermented water, organic acids, fatty acids, general oils, biodiesel for FID and other compounds containing the nitrogen groups for TEA (such as nitrosamines) etc.

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Description : It used for analyzing of volatile organic compounds, VOCs and semi-volatile organic compounds, SVOCs based on the principle of separation between two phases, consisting of stationary phase and mobile phase. When the mixtures were injected into the inlet area, which is the area that provides heat is high enough to make the liquid phase change to gas phase and the gas phase of the mixtures will move on into the column by using the carrier gas and the heating of the column with the oven, causing the separation process. Then, the mixture can be separated within the column will be entered into the detector and the result will be displayed in the form of chromatograms.

Sample type : Which can be tested both qualitative and quantitative analysis in many types of sample such as alcohols, various fermented water, organic acids, fatty acids, general oils and biodiesel, etc.

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Description: used for analyzing of volatile, VOCs and semi-volatile organic compounds, SVOCs based on the principle of separation between two phases, consisting of stationary phase and mobile phase. When the mixtures were injected into the inlet area, which is the area that provides heat is high enough to make the liquid phase change to gas phase and the gas phase of the mixtures will move on into the column by using the carrier gas and the heating of the column with the oven, causing the separation process. Then, the mixture can be separated within the column will be entered into the detector and the result will be displayed in the form of chromatograms.

Sample type : Which can be tested both qualitative and quantitative analysis in biogas samples such as hydrogen gas, nitrogen gas, methane gas and carbon dioxide gas for TCD hydrogen sulfide and other compounds containing phosphorus groups for FPD etc.

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Description: It used for analyzing of volatile, VOCs and semi-volatile organic compounds, SVOCs based on the principle of separation between two phases, consisting of stationary phase and mobile phase. When the mixtures were injected into the inlet area, which is the area that provides heat is high enough to make the liquid phase change to gas phase and the gas phase of the mixtures will move on into the column by using the carrier gas and the heating of the column with the oven, causing the separation process. Then, the mixture can be separated within the column will be entered into the detector and the result will be displayed in the form of chromatograms.

Sample type : alcohols, various fermented water, organic acids, fatty acids, general oils and biodiesel, etc.

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Description: It used for analyzing of volatile, VOCs and semi-volatile organic compounds, SVOCs based on the principle of separation between two phases, consisting of stationary phase and mobile phase. When the mixtures were injected into the inlet area, which is the area that provides heat is high enough to make the liquid phase change to gas phase and the gas phase of the mixtures will move on into the column by using the carrier gas and the heating of the column with the oven, causing the separation process. Then, the mixture can be separated within the column will be entered into the detector and the result will be displayed in the form of chromatograms.

Sample type : alcohols, various fermented water, organic acids, fatty acids, general oils, biodiesel for FID, Biogas samples such as hydrogen gas, nitrogen gas, methane gas and carbon dioxide gas for TCD hydrogen sulfide and other compounds containing phosphorus groups for FPD, etc.

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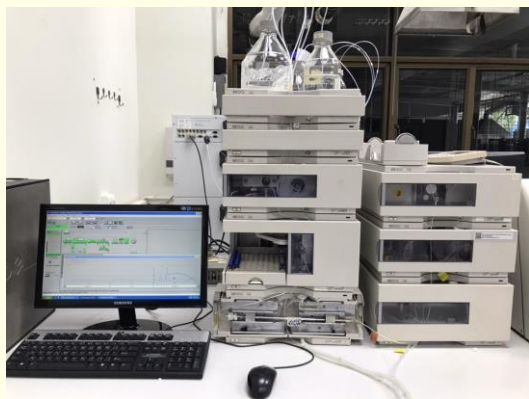
Description: Gel permeation chromatography (GPC) or Size Exclusion Chromatography (SEC) is a technique for separating the substances based on the size. The large size of polymers will move from column to the detector first due to, it cannot be facilitated into the porous of stationary phase, while the small size of polymers move from column to the detector later due to, it can be facilitated into porous of stationary phase causing the separation process within the column. This technique is often used for the analysis of molecular weight in various polymers, which can be described by a variety of definitions for molecular weight including; the number average molecular weight (M_n), the weight average molecular weight (M_w), the size average molecular weight (M_z), the viscosity molecular weight (M_v) and the polydispersity index (PDI) of molecular weight of polymers etc.

Sample type : both of natural polymers and synthetic polymers, etc.

HIGH PERFORMANCE LIQUID CHROMATOGRAPH (HPLC)-DIODE ARRAY DETECTOR (DAD), REFLECTIVE INDEX DETECTOR (RID), FLUORESCENCE DETECTOR (FLD)

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Description: HPLC works following the basic principle of thin layer chromatography or column chromatography, where it has a stationary phase and a mobile phase. The mobile phase flows through the stationary phase and carries the components of the mixture with different components travel at different rates. Thus the components separated and found in different region in chromatography to separate and detected signal with detector then identify and quantify.

Sample type: It can analyze both qualitatively and quantitatively of samples such as sugar (glucose, fructose, sucrose and xylose), preservatives, and extracts, etc.

HIGH PERFORMANCE LIQUID CHROMATOGRAPH (HPLC)-DIODE ARRAY DETECTOR (DAD), ELECTROCHEMICAL DETECTOR (EDC), EVAPORATIVE LIGHT SCATTERING DETECTOR (ELSD)

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ตัวตรวจวัด / Detector : Diode Array Detector (DAD), Electrochemical Detector (ECD),
Evaporative Light Scattering Detector (ELSD)

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Description: HPLC works following the basic principle of thin layer chromatography or column chromatography, where it has a stationary phase and a mobile phase. The mobile phase flows through the stationary phase and carries the components of the mixture with different components travel at different rates. Thus the components separated and found in different region in chromatography to separate and detected signal with detector then identify and quantify.

Sample types : It can analyze both qualitatively and quantitatively, with samples such as sugar (glucose, fructose, sucrose and xylose), preservatives, and extracts *etc.*

HIGH PERFORMANCE LIQUID CHROMATOGRAPH (HPLC)- VARIABLE WAVELENGTH DETECTOR (VWD)

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Description: HPLC works following the basic principle of thin layer chromatography or column chromatography, where it has a stationary phase and a mobile phase. The mobile phase flows through the stationary phase and carries the components of the mixture with different components travel at different rates. Thus the components separated and found in different region in chromatography to separate and detected signal with Detector then identify and quantify.

Sample types: Liquid samples such as preservatives in foods and active agents in cordyceps, etc.

HIGH PERFORMANCE LIQUID CHROMATOGRAPH (HPLC)-UV- Vis DETECTOR (UV-Vis), REFLECTIVE INDEX DETECTOR (RID)

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ตัวตรวจวัด / detector : UV-Vis Detector (UV-Vis), Reflective Index Detector (RID)

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Description: HPLC works following the basic principle of thin layer chromatography or column chromatography, where it has a stationary phase and a mobile phase. The mobile phase flows through the stationary phase and carries the components of the mixture with different components travel at different rates. Thus, the components separated and found in different region in chromatography to separate and detected signal with Detector then identify and quantify.

Sample types: Liquid samples such as preservatives in foods and active agents in cordyceps, etc.

HIGH PERFORMANCE LIQUID CHROMATOGRAPH (HPLC)-DIODE ARRAY DETECTOR (DAD), REFLECTIVE INDEX DETECTOR (RID), FLUORESCENCE DETECTOR (FLD)

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ตัวตรวจวัด / Detector : Diode Array Detector (DAD), Reflective Index Detector (RID),
Fluorescence Detector (FLD)

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Description: HPLC works following the basic principle of thin layer chromatography or column chromatography, where it has a stationary phase and a mobile phase. The mobile phase flows through the stationary phase and carries the components of the mixture with different components travel at different rates. Thus, the components separated and found in different region in chromatography to separate and detected signal with detector then identify and quantify.

Sample type: Liquid samples such as preservatives in foods and active agents in cordyceps, etc.

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Description: Fourier Transform Infrared Spectrometer use to analyze the functional group of the chemical compounds, by using the principle of absorbing the middle infrared region (400 cm^{-1} - $4,000\text{ cm}^{-1}$). When the molecule receives energy from the infrared radiation that has the frequency corresponding to the frequency of vibration or rotation of the covalent bond in the molecule, the molecule will absorb the IR radiation and change the dipole moment of the molecule. The device measures the intensity per frequency or wavelength to get the

Interferogram spectrum. This analysis results are accurate and high sensitivity with less time to analyze. This machine can compare the similarities or differences of the desired spectrum and compare the spectrum with the existing library in the device. It can measure in 2 ways: absorbance, a measurement of the samples light absorption, and % transmittance, a measurement of the light shining through the sample. The analysis techniques that are choosing depend on the property of the substances. The techniques used in the analysis are as following

- KBr Pellet
- Neat
- ATR
- Diffuse
- Photoacoustic

Sample type: The samples can be powders, solids and liquids such as synthetic substances, natural product extracts (plant, herb, animal), natural rubber, synthetic rubber, plastic and polymers, *etc.*

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Description: Fourier Transform Infrared spectrometer use to analyze the functional groups of the chemical compounds, by using the principle of absorbing the middle infrared region (400 cm^{-1} - $4,000\text{ cm}^{-1}$). When the molecule receives energy from the infrared radiation that has the frequency corresponding to the frequency of vibration or rotation of the covalent bond in the molecule, the molecule will absorb the IR radiation and change the dipole moment of the molecule. The device measures the intensity per frequency or wavelength to get the Interferogram spectrum. This analysis results are accurate and high sensitivity with less time to analyze. This machine can compare the similarities or differences of the desired spectrum and

compare the spectrum with the existing library in the device. For FTIR-Microscope can be measured both transmission and reflection modes. It can measured the very small samples.

Sample types: The samples can be powders, solids and liquids such as synthetic substances, natural product extracts (plant, herb and animal), natural rubbers, synthetic rubbers, plastics, polymers, *etc.*

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Description: Raman spectrometer is used to study the molecular structure and to analyze the functional groups. The samples are analyzed under an optical microscope that can analyze in the particle size area ranging from smaller than microns to millimeters. The spectrum analysis can be done by the laser irradiation and measuring the light scattering of the sample from the raman phenomenon. The raman spectrum is unique to the molecules and the crystal structures which can be used to identify the molecular types and evaluate the crystal structures. The machine has 3 standard lenses, including 10x, 50X and 100X.

Sample types: Powders solids and liquids such as gems, coating materials, and drugs, etc.

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Description: It is used for phase identification of a crystallinity of materials. X-ray diffraction technique is a non-destructive testing method.

Sample types: The samples are solids or viscous fluids such as powder and sheet forms, for example, clay, sand, rock forming minerals, metals, industrial materials, synthetic materials and ceramics, *etc.*

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Description: It is used for analyzing the organic chemical structures that have H and C as a main structure. The technique involves measuring the different energy levels of nuclei under the influence of magnetic fields with multinuclear broadband observe multinuclear prodigy cryoprobe and high S/N digital tuning with 1D-NMR and 2D-NMR applications.

Sample types: The samples are solids and liquids such as extracts, drugs and rubber etc. The sample must be able to dissolve in the solvent.

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Description: Bomb calorimeter is used to measure thermal energy of materials according to standard method, by burning materials with oxygen to give out heat which will be received by water tank and the increased water temperature is measured to calculate the thermal energy of materials.

Sample types: The samples can be both solids and liquids such as rice husk, oil, palm shell, coal and biomass products, etc.

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Description: CHNS/O Analyzer is used for analyzing the elements C H N S and O in samples for both liquids and solids by the method of burning Samples at high temperature into mixture gas. Then enter the column to separate and measure the amount of elements in the sample.

Sample types: The samples must be homogenous. The samples for analysis are such as chemicals, organic substance, inorganic substance, rubber, polymer, and sediment, etc.

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Description: CHNS/O Analyzer is used for analyzing the elements C H N S and O in samples for both liquids and solids by the method of burning samples at high temperature into mixture gas. Then enter the column to separate and measure the amount of elements in the sample.

Sample types: The samples must be homogenous. The samples for analysis are such as chemicals, organic substance, inorganic substance, rubber, polymer, and sediment, etc.

AUTOMATIC DIRECT MERCURY ANALYZER (DIRECT MERCURY1)

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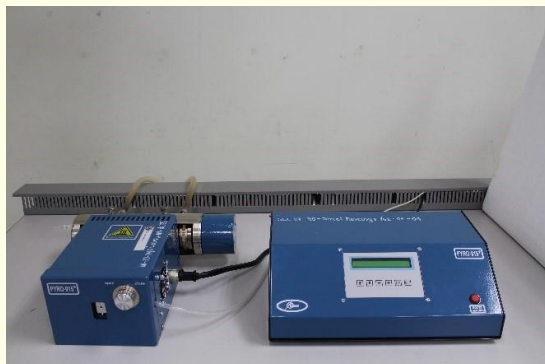
Description : It use to analyze the mercury in solid, liquid and semi-solid samples directly based on the principle of thermal decomposition by allowing mercury atoms to form free atoms to accumulate and capture gold particles into the gold amalgamation when the gold amalgamation receives heat appropriate mercury in the free-form atom moves into the sensor with the atomic absorption spectrometer (AAS) technique absorbance at 253.7 nm without sample preparation process with a measurement range of 0.001 ngHg - 2,000 ngHg.

Sample types: The samples can be both solids and liquids such as meat, soil, fertilizer, plant, food, cosmetics, plastic, medicine, oil, color, blood, urine, and waste water with homogeneous samples that contains at least 10 grams.

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Description: The Instrumentation principle is differential Zeeman atomic absorption spectrometry with High Frequency Light Polarization (ZAAS-HFM) by measurement absorption of mercury atom. The data processing system can control by using computer or standalone (in air mode).

Sample types: The samples can be both solids and liquids such as meat, soil, fertilizer, plants, foods, cosmetics, plastic, medicine, water, sea water and waste water, etc.

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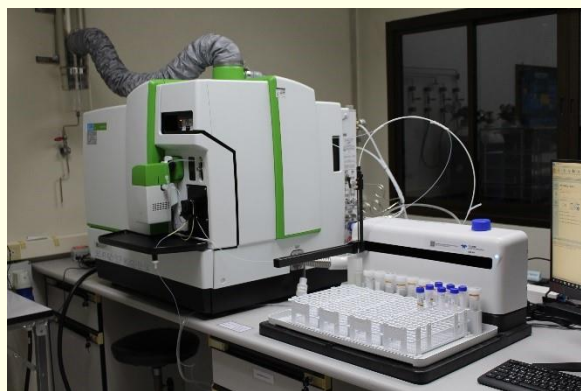
Description: Determination of mercury with the mercury system-atomic absorption spectroscopy technique (FIMS-AAS) is a technique that can measure liquid samples. Peristaltic pumps delivery the carrier and reductant. When the exact volume sample flows and react with the reductant, the reaction is mixed to be stored at the gas / liquid separator and argon gas carry $Hg^0(g)$ through the polytetrafluoroethylene (PTFE) filter to detect with sensor at absorbance 253.7 nm.

Sample type : The samples can be both solids and liquids such as meat, soil, fertilizer, plant, food, cosmetics, plastic, medicine, waste water, and water, etc.

INDUCTIVELY COUPLED PLASMA-MASS SPECTROMETER (ICP-MS)

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Description: Inductively Coupled Plasma-Mass Spectrometer (ICP-MS) is a technique to analyze the amount of elements and very small amounts of metal in ppb and ppt levels and can be analyzed in many elements at the same time by using the principle of combustion. Samples are sprayed in a spray of argon plasma to break down the charge ions + 1 and then forward to separate and measure the mass according to the mass per charge (m/z) with Mass Quadrupole type spectrometer.

Sample types: The ICP-MS can tolerate samples with a total of Total Dissolve Solid (TDS) values in approximately 1%. Various samples can be measured such as drinking water, waste water, industrial wastewater, food, plants or soil that prepared to test with controlled amount of low acid and TDS.

INDUCTIVELY COUPLED PLASMA OPTICAL EMISSION SPECTROMETER (ICP-OES1)

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Description: ICP-OES determine the concentration of elements in the ppm level.

Sample types: It can analyze both Liquid and solid samples such as fertilizer, soil, food, wood residues, rubber, and wastewater, *etc.*

INDUCTIVELY COUPLED PLASMA OPTICAL EMISSION SPECTROMETER (ICP-OES2)

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Description : MacroCHNS Analyzer is a device that analyzes the amount of carbon, hydrogen, nitrogen, and sulfur based on the principle of combustion the sample with high heat to allow the substance to be in a gas state and measure the amount of elements with the thermal conductivity detector and IR detector.

Sample types: Foods, animal feeds, dairy products, soils, plants, charcoal, and oil, etc.

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Description: OES is used to analyze and detect the amount of elements by comparison with standard substance which can be measured from ppm level to percent.

Sample types: Solid materials that are smooth surface and without pore such as steel, aluminium, copper and brass, etc.

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Description: Sulfur in Oil Analyzer is used for determination of sulfur in petroleum products. It can suggest quantity by comparing with standard substance which can be measured from the percent or ppm levels.

Sample types: Petroleum products such as biodiesel oil (B5), diesel oil, crude oil and light oil, etc.

TOTAL ORGANIC CARBON ANALYZER (TOC) for Liquid

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Description : TOC is used to analyze total organic carbon and nitrogen (TN) in the samples by combustion with high temperature.

Sample type : Liquid samples such as natural water, drinking water, and wastewater *etc.*

TOTAL ORGANIC CARBON ANALYZER (TOC) for solid

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Description: TOC is used to analyze total organic carbon, total inorganic carbon and total carbon by combustion with high temperature.

Sample types: Homogeneous solid samples such as soil, and chemicals, etc.

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Description: XRF is used to analyze elements in semi quantitation from Be to U. It can suggest quantity by comparing with standard substance which can be measured from the ppm to percent levels.

Sample types: The samples are both solids and liquids such as rock, clay, mineral, metal, and oil, etc.

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Description: Nitrogen or protein content analysis by Kjeldahl technique is performed by protein degradation with sulfuric acid. The nitrogen in the sample will be changed to ammonium sulfate $((\text{NH}_4)_2\text{SO}_4)$ under high temperature conditions with catalysts such as CuSO_4 , etc. After that ammonia is distilled out of the digested solution and is analyzed by means of titration techniques which determine the amount of ammonia in the form of ammonia compounds.

Sample types: Liquids and solids such as raw water, wastewater, feed, feed ingredients, forage, pet food, grain, cereal, meat, meat products, milk and dairy products, etc.

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Description: Nitrogen content analysis by Kjeldahl technique is performed by protein degradation with sulfuric acid. The nitrogen in the Sample will be changed to ammonium sulfate $((\text{NH}_4)_2\text{SO}_4)$ under high temperature conditions with catalysts such as CuSO_4 , etc. After that ammonia is distilled out of the digested solution and is analyzed by means of titration techniques which determine the amount of ammonia in the form of ammonia compounds.

Sample types: Liquid or solution samples such as raw water and wastewater, etc.

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Description: Coulometric Karl Fischer Titration used for measure the water content based on the principle of titration of samples containing water with iodine solution in methanol, sulfur dioxide and pyridine. Which iodine reacts with water, when the reaction completely, solution change to the brown color according to free iodine molecule. Then it measures the volume of iodine used in the titration. This apparatus is suitable for testing of water content in the range of 0.001-10 percent by weight.

Sample types: Many types of sample can be tested such as solvent, wood vinegar, oil, biodiesels and petroleum products, etc.

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Description: Volumetric Karl Fischer Titration used for measure the water content based on the principle of titration of samples containing water with iodine solution in methanol, sulfur dioxide and pyridine. Iodine reacts with water, when the reaction completely, solution change to the brown color according to free iodine molecule. Then, It measure the volume of iodine used in the titration. This apparatus is suitable for testing of water content in the range of 0.1-100 percent by weight. For the samples is cannot dissolved in solvents, such as herb powder, resin or plastic, using a device called drying oven It heats the water in the sample to steam and steam is sent to the tube for analysis in the vessel of the Karl Fischer machine. It can set the temperature from 50 °C -300 °C.

Sample types: solvent, wood vinegar, oil, biodiesels, petroleum products chemical and plastic, etc.

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Description: Spectroquant® Pharo300 can analyze the amount of substances in liquid samples such as chloride, nitrate, fluoride, sulfate, color, and turbidity, etc. The principle based on light absorption with xenon flash lamp light source generates wavelength of 190 nm – 1100 nm.

Sample types: Liquid samples such as natural water, drinking water, wastewater, and sea water etc. In the case of solid samples must be prepared before testing.

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Description : Spectroquant® Pharo100 can analyze the amount of substances in liquid samples such as chloride, nitrate, fluoride, sulfate, color, and turbidity, etc. The principle based on light absorption with halogen flash lamp light source generates wavelength of 320 nm – 1100 nm.

Sample types: Liquid samples such as natural water, drinking water, waste water, sea water etc. In the case of solid Samples, must be prepared before the test.

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Description : UV-Vis is used to measure both qualitatively and quantitatively to determine the amount of the sample and intensity by using the UV and the visible lights that pass through or absorbed. The wavelength is related to the amount and type of the sample, which is mostly organic compounds, complex compounds and inorganic compounds that can absorb in these wavelengths. It can measure according to standard methods in 3 types Abs,% T and % R. Measurement range - 3.00 - 3.00 Abs and wavelength range 190 nm - 2700 nm.

Sample types: The samples can be solids, films and liquids such as the surface of the sample, film, biocalcium, phenol, etc. The sample must be homogeneous and use a small amount.

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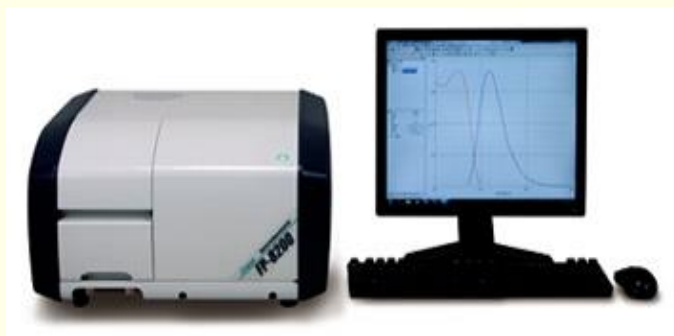
Description: UV-Vis is used to measure both qualitatively and quantitatively to determine the amount of the sample and intensity by using the UV and the visible light that pass through or absorbed. It detects light absorption value on the wavelength of 190 nm - 1100 nm. It can measure according to the standard methods in 3 types: Abs, %T and % R.

Sample types: The samples should be soluble in the organic solvent such as curcumin solution, DNA and protein, etc. It uses a small amount of the sample.

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Description: Fluorometer is used for analyzing the property of substance such as analyzing the value of excitation, emission, constant of wavelength synchronous and constant of synchronous spectral which can scan excitation and emission in 3 dimension by the method of UV absorption, causing molecule to be stimulated and shake within molecule from ground state to excited state, resulting in spectrum in the fluorescent period at the stimulated energy value of each substance.

Sample types: The samples should be solution such as toxicological study, medical study, the study of the reaction between enzymes and precursor, vitamin, and organic food substance, etc.

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Description: Titrate 794 Basic Titrino use for a testing total acid number refer to ASTM D664, total base number refer to ASTM D4739, % glycerol refer to ASTM D7637 in general oil and biodiesel. It can find the volume of titrant for titration with the sample to the end point based on the principle of potentiometric titration. It is the measurement of the electric potential that is changed while adding the volume of the titrant into the samples under continuous stirring with magnetic stirrer. The electric potential is measured as a result of the reaction between the sample and the titrant.

Sample type : General oils such as biodiesel, etc.

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Description: Total fat analyzer is used to analyze crude fat and total fats.

Sample types: The samples can be solids, liquids and semi-solids such as waste water, raw materials, intermediates and finished products in food, animal feed, and pet food, *etc.*

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Description: The analyzer can determine the protein content in food samples such as animal feed, meat, milk by measuring the only true protein and not include non-protein nitrogen. Able to determine the protein content at the range from 0.01% - 99.99% with 0.01% resolution for liquid, solid or condensed sample. It is fast analyzing process. Analyzing techniques is green technology without the use of chemicals and experimental conditions that are harmful to users.

Sample types: Liquid, solid or condensed food samples such as animal feed, meat, and milk, etc.

OXIDATION STABILITY TESTER ACCORDING TO EN 15751:2009

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Description: For Oxidation stability at 110 °C in biodiesel according to EN 15751:2009.

Sample types: General oils such as biodiesel, etc.

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Description: Fat extractor use for fat extraction by using the principle of solvent extraction.

Sample types: Solid Samples such as palm kernel, palm waste, animal feeds, etc.

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Description: Moisture analyzer can be use for determination of moisture in various types of sample. Its specifications are maximum weight capacity: 51 g and weight resolution: 0.001 g.

Sample types: Homogeneous solid samples such as chemicals, milk powder and feeding stuffs, etc.

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Description: Fiber analyzer, by digestion of sample with both acid and alkali based on AOAC method.

Sample types: Solid samples such as herbs, feeding stuffs and instant foods, etc.

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Description: Automatic BOD and PC-BOD: BOD Analyzer performs the determination of residue oxygen in water sample. The system must follow standard methods 5210 B 5 day BOD Test.

Sample types: Water such as raw water, and wastewater, etc.

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FIELD EMISSION SCANNING ELECTRON MICROSCOPE

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Description: Field emission scanning electron microscope (FESEM) provides topographical information. The highest magnification is 1,000,000x. Attach with Energy Dispersive X-Ray Spectrometer; EDS, (Oxford). EDS Analysis provides elemental analysis of a sample.

Sample types: Magnetic sample, Material sample (metal, concrete, film, rubber, and polymer, etc.), and biological sample (fungi, bacteria, pollen, and animal tissue, etc.)

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Description: The Leica M205 FCA fluorescence stereo microscope unites high-end imaging capabilities magnification 3.9-160. Besides it com imaging with fluorescence such as ETDAPT filter for exciting of UV light, ETGFP filter for exciting of blue light and ETTXR filter for exciting of greenlight.

Sample type: Biological samples: ant, polychaeta, and pollen etc., and material samples: rubber, and polymer, etc.

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Description: Scanning electron microscope is a technique for study microstructure of the sample at maximum magnification is 20,000 times. Application of SEM is for surface topography; size of particle, porous, composite materials and grain metal for material samples include muscle, fungi fiber, size and surface of pollen for biological samples.

Sample type: dry solid sample, since the specimen chamber is at high vacuum. Hard, dry materials such as wood, bone, feathers, dried insects, or shells (including egg shells) can be examined with little further treatment, but living cells and tissues and whole, soft-bodied organisms require chemical fixation to preserve and stabilize their structure.

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Description: The FEI Quanta 400 SEM is a tungsten filament SEM, highest magnification is 300,000. It can produce enlarged images of a variety of specimen achieving $\times$. The High Voltage can be changed to span any values from 200V to 30kV. Attach with Energy Dispersive X-Ray Spectrometer; EDS, (Oxford). EDS Analysis provides elemental analysis of sample.

Sample type : The samples can be material sample such as rice husk, oil, palm shell, coal or biomass product, metal, concrete, film, rubber, and polymer, etc. Biological sample such as fungi, bacteria, pollen, and animal tissue, etc.

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Description: Scanning Electron Microscope (SEM) with a tungsten electron source has both SEI (Secondary Electron Image) and BEI (Backscatter Electron Image) image detector. SEM is often used in the structural and surface analysis of biological and materials samples. The maximum sample size is 300 mm in diameter and 5 kg in weight. Including, the elemental composition analysis of sample was analyzed by EDS (Energy Dispersive X-ray Spectroscopy) technique from boron (B) to californium (Cf). The highest magnification of SEM is 300,000 times. The sample can observe in high-vacuum or low-vacuum conditions. In addition, UVD-STEM (Ultra Variable-

Pressure Detector-Scanning Transmission Electron Microscopy) technique is utilized to image the internal structure of the sample.

Sample type : Material sample (metal, concrete, film, glove, fabric, natural fiber, rubber, polymer etc.) and biological sample (Fungi, Bacteria, Pollen, Teeth, Animal tissue etc.).

TRANSMISSION ELECTRON MICROSCOPE (TEM)

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Description: Transmission electron is a microscopy technique in which a beam of electrons is transmitted through a specimen to form an image. The specimen is most often an ultrathin section less than 100 nm thick or a suspension on a grid. Transmission electron microscopes are capable of imaging at a significantly higher resolution than light microscopes, magnification up to 1,500,000 times.

Sample type: Biological samples: Plant, virus, tissue and microorganism and material samples: polymer, nanoparticle, metal and micro-emulsion.

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Description: High sensitivity XRF qualitative and quantitative analysis with an X-ray imaging microscope. Analytical procedures include multi point analysis, line analysis mapping analysis. The Sample area is a minimum of 10 micrometers by 10 micrometers and maximum Sample area of 10 centimeters.

Sample types : Solid and liquid samples such as rubber pillow, milk and plastic, etc.

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Description : AFM is a microscopy technique that can provide three-dimensional images of virtually any surface at nanometer-scale resolution. It is possible to test in both air and liquid conditions. It relies on the force between a sharp probe and the surface, which is detected while scanning the probe over the Sample. AFM can characterize a wide array of mechanical properties (e.g. adhesion, stiffness, friction, dissipation), electrical properties (e.g. capacitance, electrostatic forces, work function, electrical current), magnetic properties, and optical spectroscopic properties. ANA software (Automated Nanomechanical Analysis) can investigate the nanomechanical properties of materials such as cells, tissues, scaffolds, hydrogels, and polymers on multiple or large Samples in an easy-to-use fashion.

Sample types : Various materials e.g. cells, tissues, scaffolds, hydrogels, coating surface and polymers.

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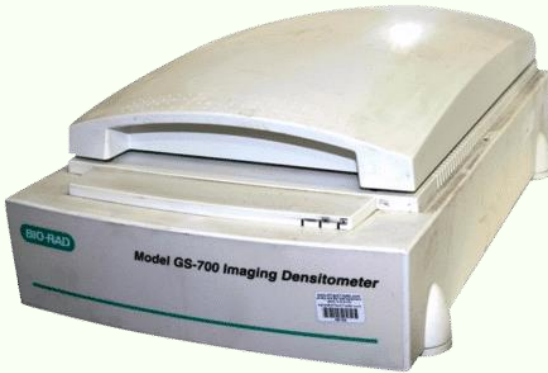
Description : The L-8900 is a high-speed amino acid analyzer with high sensitivity that is also capable of analyzing physiological fluids and protein hydrolysates. Speed Analysis Quantitative analysis of amino acids in the 30 minutes for a specify type and amount of amino acid for determine its label nutritional.

Sample types : The samples can be both solid and liquid, such as food, feed, beverage, and herbs, etc.

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Description : Densitometer can scan Electrophoresis transparent and opaque as digital images to analyze density Able to record gel film, image, analyze size, quantity and molecular mass of DNA pieces

Sample type : Magnetic sample, material sample (metal, concrete, film, Rubber, and Polymer, etc.), and biological sample (fungi, bacteria, pollen, and animal tissue, etc.)

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Description : The machine used the isolation of compounds, such as amino acids, protein ions and nucleic acids, A substance that has an electric charge moving to the polar opposite, speed of movement is depends on amount of charge on the molecule, shape and size of the molecules and electricity.

Sample type : Separate the bar, DNA, protein, PCR product

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Description : Gel documentation used to record images and positions of DNA, RNA and single stand protein from blot or gel electrophoresis. The strips of this isolated DNA must be labeled with dye. It is a fluorescence compound that glows and can absorb ultraviolet radiation and the light in the visible range to observe the appearance of the strip of DNA isolated on the gel sheet and can control the work direct computer processing

Sample types : X-ray film, autoradiograms, SDS-face gel, silver gel (wet or dry), polaroid, nitrocellulose, and ethidium bromide gels.

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Description : Instrument use to identification Microbial using Mass spectrometry and construction of protein structure. The results will be used to compare data such as library for rapid detection of bacteria.

Sample type : Food, environment, and single colony microorganisms samples.

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Description : Microplate Reader measures the absorbance of light at the end point or kinetic the reaction in the 96-Well Plate or interest wavelength. The machine can measure the absorbance at wavelengths from 200 - 999 nm.

Sample type : Juice, free radical inhibitors and substances in the immune system, *etc.*

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Description : Next generation sequencing (NGS), massively parallel or deep sequencing are related terms that describe a DNA sequencing technology which has revolutionised genomic research. To date, these technologies have been applied in a variety of contexts, including whole genome sequencing, De novo sequencing, Epigenetics, and Metagenomics to accelerate biological and biomedical research. Using the diagnosis to consider treatment.

Sample type : Analysis of plant, animals, microorganisms Samples have been extracted or DNA.

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Description : Instrument for preparing Sample molecular for amplification DNA Organisms include plants, animals, bacteria, viruses, yeast, and Recombinant DNA, *etc.*

Sample type : Analysis of plant, animals, microorganisms samples have been extracted or DNA.

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Description : Real-time PCR is a technique used in medicine as applied to detect avian flu (Influenza A H5N1) or other pathogens. Study the mutations of the gene, only one (SNP). Industry used to monitor raw materials and processed products from plants and animals that have been genetic engineering (Genetically Modified Organisms, GMOs) and measurement of DNA.

Sample type : Analysis of plant, animals, microorganisms Samples have been extracted or DNA.

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Description : A device for analyzing aflatoxin, a toxic substance produced by mycotoxin fungi based on the principle of Immunoreceptor test. The extract must be extracted from the Sample and drip onto the test kit (Strip.) Results can be quickly known in terms of quantity and quality.

Sample types : Cereal seeds, beans, dried seeds and oilseeds such as corn, sorghum, rice, wheat, peanuts, coconuts, herbs, spices, and in almost all processed products using raw materials from agricultural products contaminated with this type of fungus.

LIQUID SCINTILLATION COUNTER AND LUMINATER

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Description : Real-time PCR is technique used in medicine as applied to detect avian flu (Influenza A H5N1) or other pathogens. Study the mutations of the gene, only one (SNP). Industry used to monitor raw materials and processed products from plants and animals that have been genetic engineering (Genetically Modified Organisms, GMOs) and measurement of DNA.

Sample type : Analysis of plant, animals, microorganisms Samples have been extracted or DNA.

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Description : The Thermo Scientific NanoDrop Lite is a small, stand-alone UV spectrophotometer developed specifically for micro volume measurements of nucleotides such as nucleic acids and purified proteins. The NanoDrop™ Lite comes with preloaded software and is designed to be used without a computer.

Sample types : DNA samples extracted from plants, animals, bacteria including sample form environmental.

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Description : Multi-mode Microplate reader is a multi-mode microplate reader with advanced LVF monochromators™, highly sensitive filters, and an ultra-fast UV/VIS spectrometer. The instrument is upgradable and can be equipped with all the leading non-isotopic detection technologies: UV/VIS absorbance. Fluorescence intensity, including FRET luminescence (flash and glow), including BRET fluorescence polarization. Time-Resolved Fluorescence, including TR-FRET and alpha screen.

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Description : Compression molding is common process used for both thermoplastic and thermoset stock shape materials. Compression molding is accomplished by placing the plastic material (can be a granular or pelletized form) in a mold cavity to be formed by heat and pressure.

Sample type : Rubber compounds, plastic, composite, and thermoplastic elastomer, etc.

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Description : Compression molding is common process used for both thermoplastic and thermoset stock shape materials. Compression molding is accomplished by placing the plastic material (can be a granular or pelletized form) in a mold cavity to be formed by heat and pressure.

Sample type : Rubber compounds, Plastic, Composite, and Thermoplastic elastomer *etc.*

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Description: Internal Mixer, 120 cc, used for compounding rubber and thermoplastic elastomer. Typically rubber internal mixers consist of a completely enclosed mixing chamber in which two rotors operate. The tester can also be set rotor speed and temperature.

Sample type : Rubber compounds, plastic, composite, and thermoplastic elastomer, etc.

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Description : Internal Mixer, 300 cc, used for compounding rubber and thermoplastic elastomer. Typically rubber internal mixers consist of a completely enclosed mixing chamber in which two rotors operate. The tester can also be set rotor speed and temperature.

Sample type : Rubber compounds, plastic, composite, and thermoplastic elastomer, etc.

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Description : Rubber mixing system, open system, grinding rubber into homogeneous, mixed rubber and chemicals together.

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Description : Preparation of test pieces by cutting Sample with die cut according to ISO 37 and ASTM D412, etc.

Sample type : Vulcanized rubber, Compound rubber, Thermoplastic rubber, and plastic sheet, etc.

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Description : For preparation of rubber Samples from to thin sheets in the range of 0.5 mm - 20 mm, the surface is smooth and consistent, suitable for slides of rubber Samples with a hardness not exceeding 65 Shore A, with a resolution of 0.1 mm.

Sample type : Vulcanized rubber, various rubber sheets and other rubber products such as rubber speed glider shockproof rubber etc.

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Description : Use for preparation of cellular foam sample in any size. It has a large area to place samples (3 m²). Easy to cut sample 25 mm in wide and 50 nm in thickness. Give a smoot cutting surface and regularly. And can measure length for 0.5 mm accurated.

Sample type : Rubber cellular foam, polyurethane foam, rubber cellular foam pillow, rubber cellular foam mattress.

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Description : The latex rubber stability test machine is apparatus tests the mechanical stability of concentrated natural and synthetic latex rubber when subjected to high speed mechanical stirring or mechanical shear. Stirring is applied to the latex rubber at a constant of 14,000 rpm until the end point, or latex instability is reached. The end point is determined by the first appearance of small pieces of coagulated rubber

Sample type : Natural rubber latex and synthetic rubber latex.

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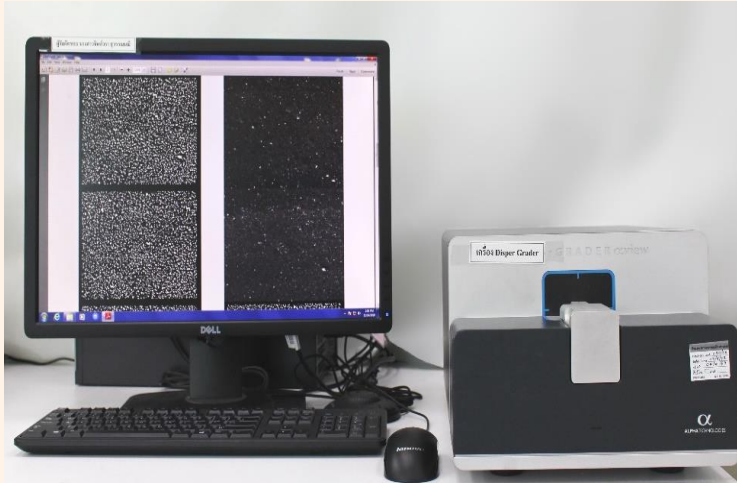
Description : A specification test for raw natural rubber and raw synthesis rubber. If keep rubber for long term they can perform oxidation reaction. Cause the Chang of Properties. The low PRI value can show low oxidation resistivity. Standard base on ASTM D3194 for this testing.

Sample type : Raw natural rubbers and raw synthesis rubbers.

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Description : Used for testing the dispersion of the mixture in the rubber compounding, capable of detecting the size, quantity, and the dispersion of composites in the compound rubber. It can analyze both white-black compound and color compound. It can also analyze the size, quantity, position of composites in the compound rubber in according to ASTM D7723.

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Description : Used for testing detecting the time when receiving the heat with rubber compound (rubber vulcanization) in according to ASTM D5289, which is important in increasing the physical properties of rubber in order to obtain the best quality rubber product. Testing temperature of RT-200 °C, strain angle: 0.2, 0.5, 1.0, 3.0 degree.

Sample type : Rubber compound (rubber vulcanization).

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Description : Used for testing the dynamic property of rubber while vulcanizing with rubber compound in accordance to ASTM D6048, ASTM D6204. It operates the duty of MDR and DMTA, capable of measuring the property of before, during, and after vulcanizing. The testing can indicate the quality control in the rubber manufacturing process in both research and industry. The frequency ranges from 0.1 to 3000, strain value $\pm 0.3\%$ to $\pm 1250\%$ (± 0.02 to ± 90 degrees), capable of detecting the G' , G'' , G^* , S' , S'' , S^*

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Description : Viscosity Test for Rubber or Rubber Compound That not vulcanized. Mostly use in industry to control rheology of rubber compound. Can Test from Room Temperature to 200 °C. The rotor rotate in 2 rpm and detect shear force from rotation resistivity of sample and rotor. Standard base on ASTM D1646 for this testing.

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Description : A capillary rheometer is an apparatus designed to measure shear viscosity and other rheological properties. Capillary rheometers for plastics are piston-die systems designed to measure viscosity of polymer melts as a function of temperature and rate of deformation.

Sample type : Plastic, composite, rubber compounds and thermoplastic elastomer, *etc.*

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Description : Use for maximum tensile strength and deformation testing of material. When apply force or deformation to test pieces and observe behavior of material. The part of machine such as pneumatic grip for holding test piece, capacity of load cell 100 N, 500 N, and 1 kN, and light extensometer to detect a deformation of each material.

Sample type : Rubber sheet, flooring rubber, rubber product, films, plastic, tire, cable, fabric, and paper. Standard for testing such as ASTM D 638, ASTM D 412, ASTM D575, ASTM D624, ISO 37, EN455-2.

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Description : Use for maximum tensile strength and deformation testing of material. When apply force or deformation to test pieces and observe behavior of material. The part of machine such as pneumatic grip for holding test piece, capacity of load cell 2.5 kN and 5 kN, light extensometer to detect a deformation of each material, and testing chamber for control temperature from -70 °C to 250 °C. Generally use for tensile strength and tear strength of plastic, film, and rubber.

Sample type : Rubber sheet, flooring rubber, rubber product, films, plastic, tire, cable, fabric, and paper. Standard for Testing such as ASTM D 638, ASTM D 412, ASTM D575, ASTM D624, ISO 37, EN455-2.

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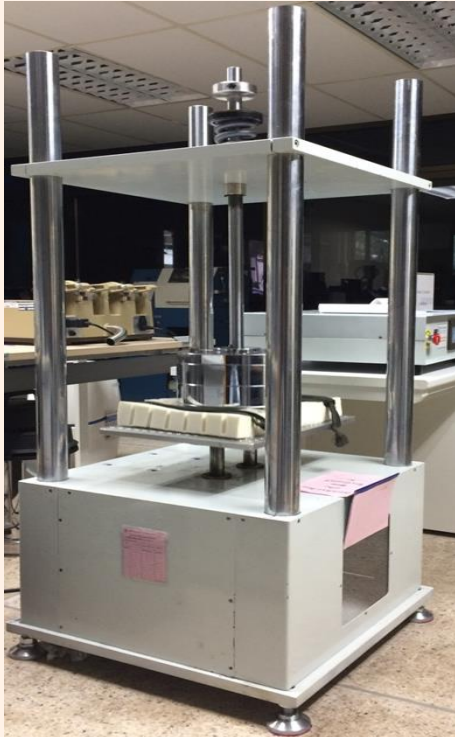
Description : For dynamic foam pounding and indentation hardness index can be before and after pounding test

Sample type : Cellular material such as latex for pillow, latex for mattress or another cellular products from polyurethane, polyethylene, synthetic fibers, glass fibers and asbestos, etc.

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Description : Foam constant load pounding machine is use to determine the thickness
reduced value and indentation force deflection decreased value of cellular material.

Sample type : Cellular material and rubber foam etc.

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Description : This universal measurement of firmness and hardness is based on a physical property called the indentation force deflection (IFD) Large Support Table allowing any test width. Conform to AS 2281, AS 2282.8, ASTM F1566, ASTM D3574 - Test B, ISO 3386, ISO 2439, BS EN 1957, GB 10807, JIS K 6400.

Sample type : Mattresses, car seats, bike seats, flexible cellular and rubber pillow, etc.

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Description : Used to test the Rebound Resilience with vulcanized rubber and various rubber products in accordance to ASTM D7121. There will be factors related to energy absorption and will be calculated from the distance of the impact and rebound. There is a ball cone of 0.2 J, hemisphere of 0.2 J, hemisphere of 0.5 J. Can be control the temperature of the specimen from room temperature to 70 °C (automatic)

Sample type : Vulcanized rubber and rubber products.

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Sample type : Shore A : Soft Rubber, Elastomer, Natural Rubber Product, Polyester, Leather, Shore D : Hard Rubber, Hard Plastic, Acrylic Sheet, Polystyrene, Shore OO : Cellular Rubber, Foam Rubber, Silicone, IRHD M (Micro IRHD) : O-Ring, Thin or Small Test Piece, IRHD N : Soft Rubber, High Elastic Material, Plastically Deformation Material, IRHD L : Cellular Rubber, Foam Rubber, Silicone, IRHD H : Hard Rubber, Hard Plastic, Acrylic Sheet, Polystyrene

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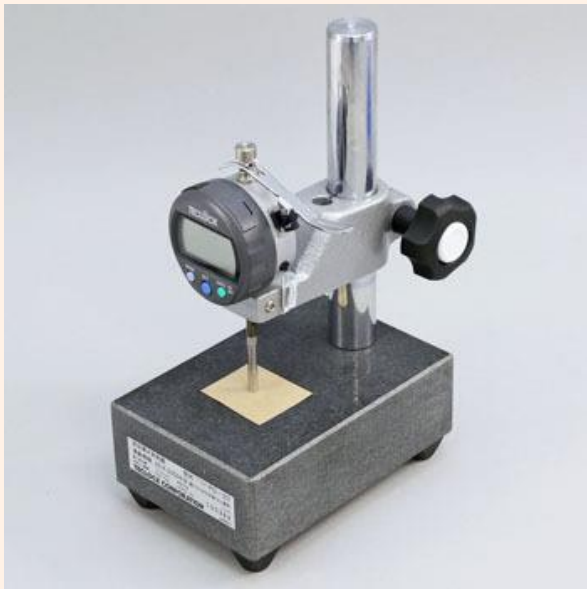
Description : Measure the *thickness* in range of 0.002 mm - 20.000 mm

Sample types : Thickness gauge is used for testing the thickness of work piece such as rubber, plastic, film, condom, and thin sheet materials, *etc.*

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Description : Measure the *thickness* in range of 0.01 mm – 10.00 mm.

Sample type : Plastic, film, rubber *etc.*

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Description : Using to Measure Density in g/cm^3 and Specific Gravity. Can Measure Accurate 0.0001 g/cm^3 . Sample size 0.001 g up to 300 and use to measure density of Rubber and Plastic.

Sample type : Rubber product and plastic.

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Description : Testing of surface friction property (Pendulum Test Value, PTV)

Sample type : Type of products such as stadium floor, multipurpose floor, rubber field, road surface, tiled floor and any pavement.

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Description : Test the air permeability in the form of air flow through the vertical Sample by measuring the air permeability can be for test between 1.4 mm / sec up to 8056 mm / sec.

Sample type : Foam, sponge and textile industry, such as woven fabric, knitted fabric, artificial leather, felt, velvet and paper *etc.*

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Description : Use for study properties of liquid on material surface. Such as Hydrophilicity Hydrophobicity, Wettability, Contact angle, Surface and interfacial tension, Surface energy of solids. Automatic Static and Dynamic Contact Angle analyzation with Sessile drop method in $0^\circ - 180^\circ$. Surface tension of liquid with Pendant drop method in range $0.01 \text{ mN/m} - 2,000 \text{ mN/m}$

Sample type : Solid material, liquid, chemical, coating material, composite material, paint, thin films, stain remover, cleanser, and cosmetic.

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Description : Low temperature compression set tester of rubber and plastic under low temperatures up to -45°C . The machine consists of 4.720 mm high spacer and can be tested according to ISO 815-2 and JIS K6261.

Sample type : Products suitable for analysis such as rubber, plastic or thermoplastic elastomers used in automotive rubber parts, etc.

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Description : Testing for property the elasticity of rubber, by finding the rubber shrinkage at low temperatures according to JIS K6261, ASTM D1329 and ISO 292. The temperature range is between -70°C to 30°C. It can test 6 work pieces simultaneously.

Sample type : Rubber tube, O-ring, tires etc.

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Description : Brittleness temperature instrument is used to measure the brittleness temperature of rubber, plastic and thermoplastic through hammering technique in the temperature range of 30°C to 100- °C following to the ASTM D746, ISO 812, ASTM D2137 and JIS K 6261.

Sample type : Rubber, plastic and thermoplastic elastomer, and automotive Rubber Parts.

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Description : Internal Mixer, 300 cc, used for compounding rubber and thermoplastic elastomer. Typically rubber internal mixers consist of a completely enclosed mixing chamber in which two rotors operate. The tester can also be set rotor speed and temperature. analysis the hardness low temperature of vulcanized rubber or thermoplastic in the temperature range of 30-70 °C in accordance with astm d1053, jis k 6261 and iso 1432, using the 80 degree twist technique. the result show the relative modulus temperature values at t2, t5, t10 and t100.

Sample type : Rubber compounds or thermoplastic used in automotive rubber parts.

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Description : Disc shaped test piece at a constant speed, to which an abrasion wheel is pressed with a fixed load. The abrasion wheel is set at an angle to the test piece and weight loss will be observed.

Sample type : Vulcanized rubber and plastic sheet coating etc.

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Description : Wear endurance tester type NBS abrasion according to ASTM D163 can test 3 work pieces simultaneously. Sample size for thickness 3.5 mm - 10.00 mm

Sample type : Rubber compound, rubber insole, rubber heel shoes, and vulcanized rubber, etc.

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Description : Taber is used for testing the wear of materials according to ASTM D4060. It consists of testing head, 500 gram and 1,000 gram and polishing stone of standard.

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Description : Used for testing the abrasion resistance of DIN with vulcanized rubber, rubber products in according to DIN 53516 and ISO 4649. Can be the weight selectable for 20 N maximum and the testing range can be set 20 m to 40 m.

Sample type : Vulcanized rubber and rubber products.

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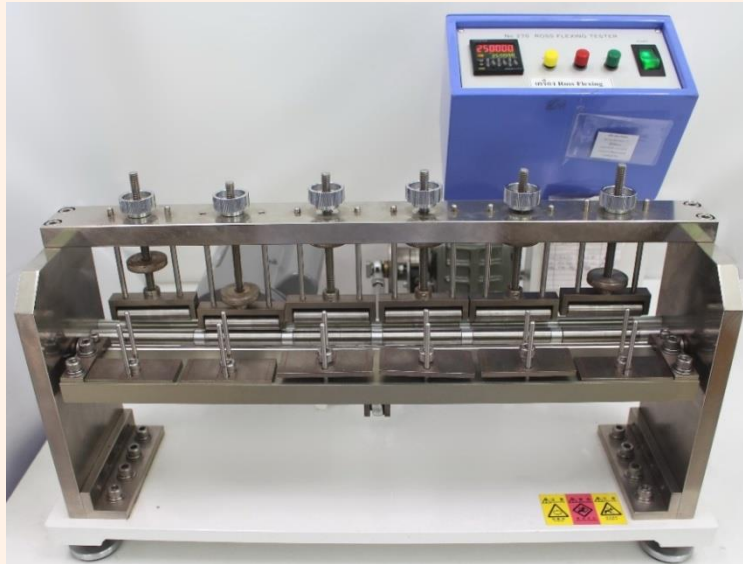
Description : Used for testing the bending in vertical repeated with speed 300 revolutions per minute is for examining the expansion of the cracks for the vulcanized rubber in according to ASTM D813.

Sample type : Vulcanized rubber and rubber products examples of types include rubber tubes, various rubber sheets.

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Description : The Ross Flexing Tester is used for testing the bending in horizontal plane is repeated in the direction of 90 degree with speed 100 revolutions per minute is for examining the expansion of the cracks for the vulcanized rubber in according to ASTM D1052. Examples of types include slippers, rubber boots, etc.

Sample type : Vulcanized rubber and rubber products examples of types include slippers, rubber boots, etc.

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Description : Compression set tester is used for measuring the compressive stresses property of materials. It can be placed in chamber with the standard Hi-Temp or Low-Temp according to ASTM D395, ISO 1856Method B, ISO 815

Sample type : Vulcanized rubber, Rubber foam *etc.*

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Description : Used to test stress relaxation with rubber and plastic in according to DIN 53504 and ISO 527. Report test results of temperature limits, TSSR-Index, and Crosslink density

Sample type : Rubber, plastic and including O-rings, rubber tubes, and valves, etc.

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Description : Tester is used to measure the increasing creep by applying static load to plastic test specimens. By choosing the Chuck, the tester enables to conduct tensile test, condensation test, and bending test. The tester can also be assorted with test functions such as relaxation test, heat shrinkage test.

Sample type : Rubber and thermoplastic elastomer *etc.*

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Description : Goodrich Flex Meter Flexometer use for measurement the heat built-up and deformation of rubber compounds. This test, which is normally performed at 30 Hz, a 1 MPa static load followed by a dynamic stroke of 4.45 mm, measures the temperature rise, creep and permanent deformation of a rubber material under dynamic load during and after the test. The test is based on the standards DIN 53 533, part 3, ISO 4666/3, (ISO 4666/4), BS 903, part A50 or ASTM D 623.

Sample type : Rubber and thermoplastic elastomer, etc.

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Description : Weathering tester expedites and controls the air condition, modeling the natural exposure for testing the quality and endurance to work piece air condition or the materials needing the endurance test such as fading, yellowing, cracking, hazing according to ASTM G155.

Sample type : Plastic bottle, Package, Film, Tile, Paint, Sanitary ware, Floor mat, Rubber block, Sports ground etc.

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Description : Accelerated weathering tester similar as weather with Xenon arc lamp. For sun light durability test and humidity durability test in dew and rain. degradation of sample shows color change, color fading, yellowing, peeling, and hazing.

Sample type : Plastic material, rubber material, woven, bottle, and packaging, etc.

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Description : Ozone Resistant Base on ASTM D1149, ISO 1431-1 Use to Measure Deteriorate of Rubber and Thermoplastic. Observe Cracking of Test Piece after Stretch in Ozone Chamber and Conditioned. The Condition Can Set to 70 °C and Ozone Concentration to 250 pphm. Deformation the piece in static and dynamic. Rotation of Holder is 2 rpm.

Sample type : Rubber product and plastic.

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Description : The resistance to multi-cell aging of vulcanized rubber and thermoplastic rubber under controlled temperature and specific time according to JIS K6257, ISO 188 and ASTM D865. It consists of 24 test tubes and can be set temperature for testing maximum to 300°C

Sample type : Vulcanized rubber, Rubber gloves, Film, Composite materials and etc.

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Description : Temperature control chamber use for high temperature resistivity of rubber and plastic. Holder rotate and control air exchange rate. Can set temperature to 300 °C and holder can rotate 5-10 rpm. Standard base on ISO 188, ASTM D573 for this testing.

Sample type : Rubber product and plastic.

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Description : Test the influence of temperature and humidity on the properties, function and lifespan of your products. Reproducible, certified and under accelerated conditions.

Sample type : Rubber product and plastic.

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Description : Can be Test Electrical resistivity up to 2×10^{16} ohms

Sample type : Rubber sheet, silicone, plastic sheet (flat surface)

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Description : It is used for testing the oxygen index for ignition with rubber, plastic and textile in according to ASTM D2863.

Sample type : Rubber, plastic and textile

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Description : Used for testing the flammability of electrical products such as wires, heat insulators, biomass materials, vulcanized rubber, plastics and polymers in accordance to ISO 5660 and ASTM E1354. Rate of heat release, Time to ignition, Mass loss rates, Smoke release rates, Thermal effects due to Effective heat of combustion, Rates of release of toxic gas such as carbon dioxide The results of such tests can be used in terms of materials research. Or use to create mathematical models to predict the severity of a burning fire

Sample type : Electrical products, biomass materials, vulcanized rubber, plastics and polymers.

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Description : Flammability tester is used to measures the fire-resistant properties of material. The test methods are based on UL 94 standard with horizontal and vertical flaming tests and associated international standards. These are 5V, VTM, HBF, HB, JIS K6911, ASTM D635, ASTM D3801, ASTM D4804, ASTM D5025, ASTM D5207 and ISO 1210.

Sample type : Rubber products and plastics.

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Description : Determines the comparative burn rates and burn resistance of textiles, plastics and other automotive interior materials. The tester conforms to Federal Motor Vehicles Safety Standard No. 302 and is also suitable for use with ISO 3795, ASTM D5132, etc.

Sample type : Products which consist of plastic, rubber and fabric including upholstery, floors, carpets, roof lining, door linings, shading etc.

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Description: Microwave digestion and extraction system is use for extraction and digestion samples before analysis by instrument such as AAS, GF-AAS, ICP, HPLC and GC. It is controlled by microprocessor system under high pressure condition by generation equal microwave energy to all tubes from magnetron with minimum frequency of 2450 MHz and maximum power is not less than 1800 watts.

Sample type: Solution sample, solid samples such as soil, wastewater, meat, and herbs, etc.

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Description: It use for drying the sample that contain water or soluble in water by using the technique that cool the sample until it freezes under vacuum condition to preserve the property of the specimen reduce the oxidation reaction and other processes that cause to destroy the sample.

Sample type: the sample can be dried in both solid and liquid, such as natural extracts, film, meat, animal feces, useful bacteria, etc.

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Description: It use for drying the sample that contain water or soluble in water. by using the technique that cool the sample until it freezes under vacuum condition to preserve the property of the specimen reduce the oxidation reaction and other processes that cause to destroy the sample.

Sample type: The sample can be dried in both solid and liquid, such as natural extracts, film, meat, animal feces, and useful bacteria, etc.

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Description : Use for drying the sample that contain water or soluble in water. by using the technique that cool the sample until it freezes under vacuum condition to preserve the property of the specimen reduce the oxidation reaction and other processes that cause to destroy the sample.

Sample type : the sample can be dried in both solid and liquid, such as natural extracts, film, meat, animal feces, and useful bacteria, etc.

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Description : The SK-300SII is a high speed auto mixer, homogenize the sample. This machine performs mixer and degassing by combining revolution and rotation movements.

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Description : Homogenizer is a multi-tasking device that can achieve high throughput sample loading task. It can automatically process up to 32 samples each run and automatically cleans the dispersing head with aqueous or solvent. The closed system reduce the noise less than 65 DB and keep you away from the solvent vapor by active exhaust system to exhaust organic volatile providing a safe environment for work.

Sample types : High humidity sample such as vegetables and fruits, high fat samples such as Meats, both cooked and raw, and high-fiber examples such as Pet food, etc.

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Description : The Waters® MV-10 ASFE® System is the first supercritical fluid extraction (SFE) system offering accelerated, multi-vessel sample extraction, resulting in a faster, greener, more selective alternative for analyze extractions from a wide variety of sample matrices.

Whether your workflow requires enriching target analyze for analytical purposes, stripping unwanted material from a product, or collecting valuable essential oils from natural products, the MV-10 ASFE System is the most efficient extraction tool available.

System features:

- Uses environmentally friendly CO₂ instead of toxic organic solvents
- Improved extraction efficiency
- Better selectivity

- Increased speed
- Higher yield
- Reduced operating costs
- ChromScope™ Software makes routine analysis, reporting, and fraction collection easy with an intuitive user-friendly interface

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Description: The 6875 Freezer/Mill is a larger cryogenic grinder with a self-contained liquid nitrogen tub and insulated case. It has one dual purpose chamber that precools and grinds samples from 0.1 to 100 grams. The grinding chamber holds one large grinding vial, one mid-sized vial, four small grinding vials, or four microvial sets.

Sample type: Bone, teeth, feeding stuffs, corn seed and plastic *etc.*

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DESCRIPTION: A sample preparation machine with automatic solid phase extraction technique is supports at least 24 samples continuous. The commonly available SPE cartridges sizes are consist of 1, 3, 6 and 12 mL. A SPE technique is a popular extraction to eliminate distractions from complicated samples to obtain substances that are of more pure interest before being measured with advanced tools such as GC, HPLC, LC / MS and GC / MS etc. This machine suitable for wide range of testing as standard methods and research.

Sample type: Sample preparation for the analysis of important substances or pesticide residues in the liquid state such as food, vegetables, fruits, drinking water, natural water, waste water and urine etc.

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290 Soi 17, Rama 9 Road, Bangkapi, Huaykwang,
Bangkok 10310

U.P.Marketing General Supply Co.,Ltd.

138, 138/11, Soi Charansanitwong 22, Charansa-
nitwong Road, Banchanglhor, Bangkok 10700

Color Global Co.,Ltd.

99/349 Moo 2, Cheangwattana Road, Tungsonghong,
Laksi, Bangkok 10210

Crest Nanosolutions (Thailand) Co., Ltd.

807/89, Phummarin Village, Moo 8, Khukot Subdis-
trict, Lam Luk Ka District, Pathum Thani 12130

Bruker Biospin AG

41 Lertpanya Building, 14Fl., Soi Lertpanya, Thanon-
Phayathai, Rajathewee, Bangkok 10400

IDI (THAILAND) COMPANY LIMITED

7/129 Central Pinklao, Office Tower A, Room 1601B
16Fl, Boromratchonni Road, Arun Amarin, Bangkok
Noi, Bangkok 10700

Chareon Tut Co., Ltd.

189 Moo 1 Theparak Road, Bangsaothong Subdistrict,
Bangsaothong District, Samutprakarn 10540

SciSpec Co., Ltd

4/11 Moo 7, Soi Bangkhae District Office, Kanchana-
pisek Rd., Bangkhae, Bangkok 10160

Miningthai Industry Council

222/2 Soi University of the Thai Chamber of
Commerce, Vibhavadi-Rangsit Road, Huai Khwang
Subdistrict, Huai Khwang District, Bangkok 10400

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- Polymer, Rubber, and Food Processing Machinery
- Block Rubber Testing Equipment